

SmartMicroScope iGO Quick Start Guide

iTunes Install:

1. Visit the iTunes Store and search for the APP “**Wi-Viewer**”
2. Install “**Wi-Viewer**”


Android Install:

1. Visit the Google Play Market and search for the APP “**Wi-Viewer**”
2. Install “**Wi-Viewer**”

NOTE: Please charge batteries for eight hours before the first use.

- A. Open the battery door on the SmartMicroScope iGO and make note of the SSID.**
- B. Turn the SmartMicroScope iGO on and leave on for about 30 seconds.**
- C. Turn the tablet on and go to SETTINGS and choose WiFi.**
- D. Choose the network SSID from the inside battery cover.**
- E. Password is “12345678”**
- F. After connected, go to the HOME screen and open the application.**

***** iOS devices ONLY: Touch the ON button on the bottom left of the screen to get the live image.

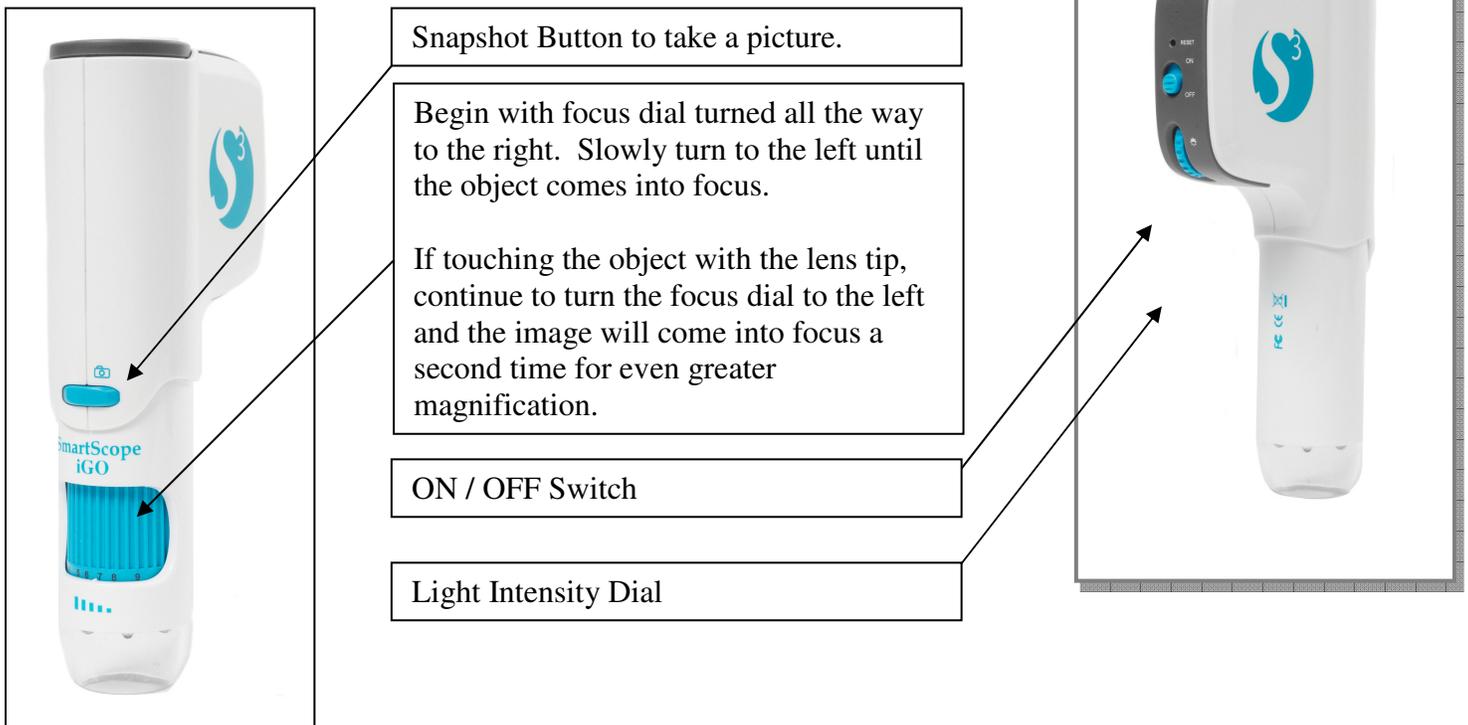

*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com